2025/05/18 11:48 1/2

Super-LOTIS Camera Pump-Down Procedure

The Vacuum pump and accessories are stored on a cart in the warm room. The pumping system consists of an Alcatel Drytel 31 pump, stainless steel vacuum hose, MKS 943 cold cathode pressure controller and a cold cathode gauge. There is also a power supply cable used to open the camera vacuum valve.

- 1. Turn telescope power on, and move the scope to El 90, Az 180 (zenith). Turn telescope power and camera power off.
- 2. Put the vacuum pump cart in the dome so the vacuum hose can reach the CCD camera.
- 3. Move the black shroud out of the way to access the bottom of the CCD camera.
- 4. Use an extension cord and a power strip to plug the vacuum pump and the gauge controller (MKS 943 cold cathode pressure) into a UPS power outlet.

On the bottom of the camera are two connections: The vacuum connection (labeled "VAC") is a smooth 1/4" diameter tube and the vacuum valve control (labeled "VAC Valve") is a phono-style power plug.

- 1. Install the compression connection on the end of the vacuum hose to the VAC inlet, and tighten.
- 2. Turn on the vacuum pump. There is a single power switch on the front, all operations are automatic. When the pump is up to speed, the green light will turn on.
- 3. After the pump is up to speed, turn on the gauge controller (black power switch on MKS 943) and then flip the small silver switch to "HV On". The digital readout will show the vacuum in a few seconds.
- 4. If the vacuum is not already at or below the 10E-3 range, then you have a leak, probably at the camera connection. Turn off the HV on the gauge controller, then turn off the gauge controller power, then turn off the vacuum pump. Reseat the vacuum connection and go back to step 8.
- 5. If the vacuum is dropping into the 10E-4 range, then turn off the HV to the gauge controller.
- 6. Plug the power cable into the VAC Valve on the camera, and use the screw threads to tighten it snugly. Plug the power supply into the power strip. THIS OPENS THE INTERNAL VACUUM VALVE IN THE CAMERA, AND EXPOSES IT TO VACUUM.
- 7. Wait 20 seconds or so, and turn on the gauge controller (HV On). It should be reading low 10E-3 and dropping.
- 8. Let the system pump for at least 1 hour.

After 1 hour or more:

- 1. Unplug the VAC Valve cable from the power strip. Unplug it from the back of the camera.
- 2. Turn the HV Off and turn the gauge controller off. Turn the pump off.
- 3. Wait for the pump to spin down. Unplug the two power cables from the power strip.
- 4. Disconnect the vacuum line from the back of the camera.
- 5. Place the pump and accessories back in the warm room.
- 6. Replace the shroud around the CCD camera.
- 7. Turn the mount power on, and the camera power back to Remote. Stow the telescope. Turn the mount power to Remote. Turn off lights and close doors.

Last update: 2019/02/26 public:kitt_peak:superlotis:superlotis_pumpdown_procedure https://lavinia.as.arizona.edu/~tscopewiki/doku.php?id=public:kitt_peak:superlotis:superlotis_pumpdown_procedure&rev=1551198725 09:32

From: https://lavinia.as.arizona.edu/~tscopewiki/ - **MOON**

Permanent link: https://lavinia.as.arizona.edu/~tscopewiki/doku.php?id=public:kitt_peak:superlotis:superlotis_pumpdown_procedure&rev=1551198725

Last update: 2019/02/26 09:32